

ILLUM-A-FIELD™

Instruction Sheet



Charging Power Pack:

1. Charge battery unit for 12 hours prior to initial use.
2. Plug charger into 110 volt AC wall outlet.
3. Plug charger cord into charge jack.
4. Small red LED (charge indicator) will glow when unit is charging.

Operating Illum-A-Field Instrument:

1. Unplug Illum-A-Field instrument from charger.
2. Plug Illum-A-Field instrument into instrument jack and switch unit on.

Switch (charge position)

Charge Indicator

Instrument Jack

Charge Jack



Do's and Don'ts:

1. Keep unit clean and dry.
2. Instruments may be cleaned with soap and water, and/or disinfectant.
3. Do NOT autoclave any component of Illum-A-Field System.
4. These instruments are currently not designed for surgery—use in the laboratory only.
5. Do NOT use these instruments with any other battery or power pack. There are internal electronics to protect the LED. Use of any power pack other than supplied by IPC will void warranty.
6. DO NOT stare into the LED unit.
7. Power Pack will power the instrument for 24 hours or more of continuous use, charge regularly.



Distributed By:

NEWCOMER SUPPLY®
Your Histology Supplier

Middleton, WI • (800) 383 • 7799 • www.newcomersupply.com